



neolab Migge GmbH
Rischerstr. 7-9
69123 Heidelberg
Deutschland
+49 (0)6221 /
8442-44
<https://www.neolab.de>
e

Umsatzsteuer-
Identifikationsnummer
:
DE 143 450 657



neolab® AXION® EyePod eye shower for taps

€461.00
plus VAT &
Shipping

Product Images



Description

The **neoLab^{AXION®} eyePod eye shower** is the ideal solution for industry and laboratories. Wherever chemicals and other substances are worked with that pose a potential hazard to humans and the environment, eye showers are a must.

Easy mounting

The eyePod eye shower can be quickly and easily mounted to standard faucets without interfering with the operation of the faucet. Simply unscrew the top of the faucet and screw the eye shower firmly onto it.

Easy to use

When foreign matter enters the eye, the faucet must be turned on and the front part of the eyePod eye shower must be rotated 180°. The eye shower is immediately active and can be used.

Increased safety

The flow and spray technology of ^{AXION®} technology also ensures increased safety and better cleaning of the eye compared to conventional eye showers. Reverse wash streams clean the eyes by gently removing debris from inside the eye and nasal cavity.

Scald protection

A built-in thermostat ensures that the water flow is stopped immediately when the water temperature reaches 37 °C. This prevents scalding of the eyes. As soon as the temperature drops again, water can flow again.

Additional Information

No.	2-1157
Manufacturer (Brand)	neoLab
EAN	4058072155698
Transport temperature	Room temperature
Showers and eye wash properties	Water jet adjustable
Connection water	3/8"
Minimum flow pressure	1 bar
Volume flow	44654

